

Thermo Scientific PetroilQuant for ARL PERFORM'X spectrometer

X-Ray Fluorescence Analytical Package

PetroilQuant™ is a comprehensive analytical package which covers the widest range of elements and concentrations in a variety of petroleum industry products. In conjunction with the Thermo Scientific™ ARL™ PERFORM'X spectrometer, PetroilQuant offers the most cost effective solution for any laboratory dealing with petroleum products.



As quality and environmental demands on the petroleum industry become stricter, key elemental contaminants require ever-lower levels of quantification, for example, stricter regulations on sulfur (S) in vehicle fuels, reduced trace metals that poison catalysts and induce corrosion such as nickel (Ni) and vanadium (V), and reduced catalyst fines aluminum (Al) and silicon (Si). The proven ability of Wavelength Dispersive X-ray Fluorescence (WDXRF) to produce highly reliable and repeatable results is increasingly solicited for such analyses in the petroleum industry. Its main advantages for such analyses are:

- Excellent repeatability
- Excellent resolution, especially for light elements (Na to Ca)
- High dynamic range (sub-ppm to 100 %)
- Flexibility in terms of measurement of analytical lines, background positions and internal references
- Little or no sample preparation in most cases (samples can be directly measured without dilution)
- Multi-element and multi-matrix capability

Numerous ASTM, ISO, IP and other international standard test methods have

been developed for petroleum industry using WDXRF. Our company offers a range of X-ray instruments and solutions depending on the elements needed, their concentration ranges, the variety of samples (liquids and solids) and the throughput requirements. The advanced ARL PERFORM'X spectrometer (Figure 1) provides full capabilities for demanding applications such as trace contaminant quantification in heavy fuels, wear metals in used lubricants, and fresh and spent catalyst characterization. The instrument is noted for high sensitivity and stability across a wide range of petroleum applications.

To provide faster solutions and meet tougher requirements in the evolving industry, we developed PetroilQuant, a unique pre-calibration program with petroleum analysts in mind. PetroilQuant is designed for use in conjunction with the ARL PERFORM'X X-ray spectrometer and is capable of quantifying up to 30 elements in a variety of light and heavy fuels, lubricants, crude oil and other petroleum process streams both quickly and cost-effectively.



Figure 1. ARL PERFORM'X WDXRF spectrometer

Thermo
SCIENTIFIC

With PetroilQuant, Thermo Scientific X-ray spectrometers are calibrated for a series of elements in a wide dynamic range covering multiple petroleum matrices. PetroilQuant requires no analytical expertise and provides a “ready to run” analytical package for beginners as well as advanced users of our XRF instruments. Further refinement or fine tuning for specific elements can always be achieved either in the factory prior to shipment or after installation using the customer’s own set of standards.

In addition to sulfur in light and heavy fuels, PetroilQuant covers other elements needed for blending control of fresh lubricants such as Zinc (Zn), Phosphorus (P), Chlorine (Cl), Calcium (Ca), Magnesium (Mg), Molybdenum (Mo), Silicon (Si), Barium (Ba) and others. The salient features of this unique package are:

- The pre-calibration program facilitates quantitative analysis for a number of elements in automotive fuels, lubricants, heavy residual oils and wear metals in lubricants
- PetroilQuant is developed with built-in knowledge and expertise in order to help the analyst deal with a variety of petroleum samples in wide concentration ranges
- PetroilQuant provides a “ready to run” analytical package for beginners as well as advanced users of our XRF instruments
- PetroilQuant saves time and money as it offers a global calibration program which otherwise would require many standards, analytical expertise and instrument time
- In conjunction with ARL PERFORM’X Series, it offers the most cost effective solution for any laboratory dealing with petroleum products

PetroilQuant can be installed on ARL PERFORM’X Series for an extended range of elements, lower concentration limits, higher throughput and faster speed of analysis.

Specifications

A series of reference materials are used to derive calibration working curves for analysis of additives in a petroleum base. This factory pre-calibration can be optimized on-site using customer’s own standard samples.

The standard PetroilQuant package for ARL PERFORM’X spectrometer includes up to 30 elements in petroleum products.

Analysis times typically range from 8 to 38 seconds depending on the element, the precision required and the power level. Table below shows typical limits of detection which are 3 sigma values from running a repeatability test on a blank oil sample.

A working curve is established for each element using the Multi-Variable-Regression incorporated in the instrument software package. This pre-calibration includes a set of three setting-up samples for maintenance of the calibration curves over time (SUS Petro1 S703976, SUS Petro2 S703977 and SUS Petro3 A036472). No standard samples are delivered with this pre-calibration.

Pre-calibrated elements

Element	Line	Analytical Range (%)	Typical LoD (ppm)	Count times 4200W (sec)	Count times 2500W (sec)	Count times 1500W (sec)
Ag	K α	LoQ-0.06	5.8	12	17	29
Al	K α	LoQ-0.06	4.7	16	23	38
As	K β	LoQ-0.06	2.7	12	17	29
Ba	L α	LoQ-0.8	1.5	8	12	19
Bi	L α	LoQ-0.12	1.3	8	12	19
Br	K α	LoQ-0.12	0.7	8	12	19
Ca	K α	LoQ-3.0	0.7	8	12	19
Cd	K α	LoQ-0.06	7.2	12	17	29
Cl	K α	LoQ-2.0	3.2	8	12	19
Co	K α	LoQ-0.06	0.6	8	12	19
Cr	K α	LoQ-0.06	0.5	8	12	19
Cu	K α	LoQ-0.12	0.4	8	12	19
Fe	K α	LoQ-0.12	0.8	8	12	19
Hg	L α	LoQ-0.06	1.4	8	12	19
K	K α	LoQ-0.8	1.5	8	12	19
Mg	K α	LoQ-0.4	12.2	18	26	43
Mn	K α	LoQ-0.06	0.6	8	12	19
Mo	K α	LoQ-0.4	1.5	8	12	19
Na	K α	LoQ-0.4	81.7	20	29	48
Ni	K α	LoQ-0.06	0.5	8	12	19
P	K α	LoQ-0.5	1.3	12	17	29
Pb	L β	LoQ-0.12	1.3	8	12	19
S	K α	LoQ-5.0	1.3	10	14	24
Sb	L α	LoQ-0.06	1.8	12	17	29
Si	K α	LoQ-0.4	2.4	14	20	34
Sn	L α	LoQ-0.06	1.5	12	17	29
Ti	K α	LoQ-0.06	0.6	8	12	19
Tl	L α	LoQ-0.06	1.2	8	12	19
V	K α	LoQ-0.06	0.5	8	12	19
Zn	K α	LoQ-0.5	0.5	8	12	19

LoQ = 3 x LoD = Limit of Quantification
LoD = Limit of Detection

www.thermoscientific.com/performx

©2013 Thermo Fisher Scientific Inc. All rights reserved. ISO is a trademark of the International Standards Organization. All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.



Thermo Fisher Scientific (Eublens)
SARL, Switzerland is ISO certified.

Africa-Other +27 11 570 1840
Australia +61 3 9757 4300
Austria +43 1 333 50 34 0
Belgium +32 53 73 42 41
Canada +1 800 530 8447
China +86 10 8419 3588
Denmark +45 70 23 62 60
Europe-Other +43 1 333 50 34 0
Finland/Norway/Sweden
+46 8 556 468 00
France +33 1 60 92 48 00
Germany +49 6103 408 1014
India +91 22 6742 9434
Italy +39 02 950 591

Japan +81 45 453 9100
Latin America +1 561 688 8700
Middle East +43 1 333 50 34 0
Netherlands +31 76 579 55 55
New Zealand +64 9 980 6700
Russia/CIS +43 1 333 50 34 0
South Africa +27 11 570 1840

Spain +34 914 845 965
Switzerland +41 21 694 71 11
UK +44 1442 233555
USA +1 800 532 4752

PS41286_E 06/13C

Thermo
SCIENTIFIC
Part of Thermo Fisher Scientific